

30 minutes Recap : Food Ingredient Analysis by LC-MS

An Advanced Analytical Technique for Food Safety and Quality

PRESENTED BY

Pongsagon Pothavorn, Sci Spec Co., Ltd.



Primary Ingredients : The Foundation of Food



Grains & Staples

Rice, flour, and sugar form food bases in many cultures.

Proteins

Meat, fish, and eggs provide essential amino acids.

Produce

Vegetables and fruits deliver vitamins and flavor diversity.

Dairy & Legumes

Milk products, legumes, seeds, and nuts support nutrition and texture.

Food Additive : Enhancing & Preserving



Type	Examples	Common Codes
Preservatives	Sodium benzoate, Sorbic acid	E211, E200
Coloring Agents	Synthetic dyes, Beta-carotene	E102, E160a
Flavourings	Vanillin, Fruit extracts	Often not coded
Sweeteners	Aspartame, Sucralose	E951, E955
Stabilizers & Thickeners	Gelatin, Gum arabic, Modified starch	E441, E414, E1400–1450
Gelling Agents	Agar, Carrageenan	E406, E407
Acidity Regulators	Citric acid	E330

Processing Aids



Enzymes

Amylase and protease aid in food processing but do not remain in products.

Solvents

Ethanol and others help extract components during manufacturing.

Filtering Agents

Used to clarify liquids; mostly removed before final food packaging.

Seasonings and Flavorings



Sauces

Soy, oyster, and fish sauces boost umami.



Basic Seasonings

Salt and sugar balance flavor profiles.

Flavourings

Natural or artificial agents enhance taste.

Herbs and Spices

Garlic, pepper, and others add aroma and zest.

Food Allergens



- Gluten from wheat, barley, rye
- Milk and dairy
- Eggs
- Peanuts and tree nuts (almonds, walnuts)
- Soybeans
- Fish and shellfish like shrimp, crab

Functional Ingredients



Functional Ingredients

- Dietary fiber boosts digestion 
- Probiotics support gut health 
- Fortified vitamins and proteins enhance nutrition 

Why it matter



Safety & Quality Assurance

Ensures ingredients meet safety standards and quality expectations.

Product Innovation

Adjust taste, texture, shelf-life. Use of novel or substitute ingredients

Label Accuracy

Verifies nutritional content, allergens, additives claims on packaging and labels.

Understanding LC-MS



Wide Compound Range

Identifies and measures diverse chemical compounds in food.

Trace Detection

Finds contaminants and residues at extremely low levels.

Minimal Preparation

Efficient analysis with less elaborate sample processing.

Case Studies: Real-World Applications



Adulterant Detection

Melamine in milk detectable at parts-per-billion levels.

Vitamin Quantification

Accurate measurement of vitamin D in fortified foods.


Food Allergens

Detection of allergens such as proteins in tree nuts, eggs, milk, soy, gluten, etc

Fatty Acid Analysis

Profiling omega-3 and omega-6 contents in oils.

Orbitrap : High-Resolution Accurate Mass

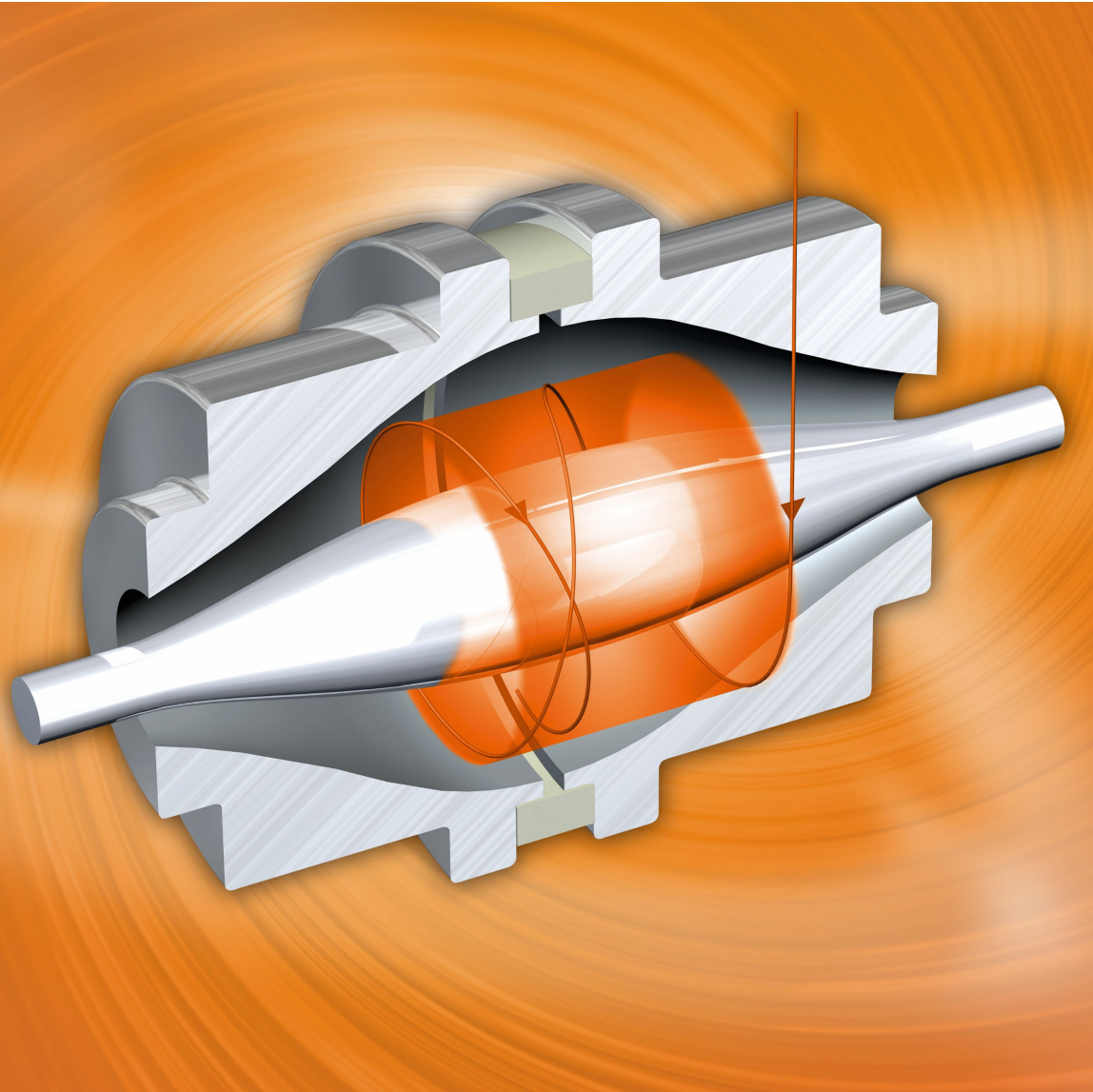

Orbitrap Analysis: Targeted & Untargeted Food Ingredient Screening

Explore how Orbitrap revolutionizes food safety and quality control. This presentation covers HRAM spectrometry principles, challenges in food matrices, and advanced screening methods.

Discovery Components



Comprehensive Profiling

Detects all compounds without prior knowledge.

Advanced Software Tools

Data analysis with XCMS and MZmine for peak detection.

I.D. & Applications

Detect adulterants, authenticate products, and assure quality.

Nutritional Analysis



Nutrient Quantification

Measures vitamins, amino acids, and essential nutrients accurately.

Label Verification

Ensures nutritional claims on packaging are truthful and reliable.


Precision

LC-MS delivers reproducible and sensitive results for nutrition.

Example

Vitamin D levels in fortified milk precisely analyzed by LC-MS.

Predicted Elemental Composition

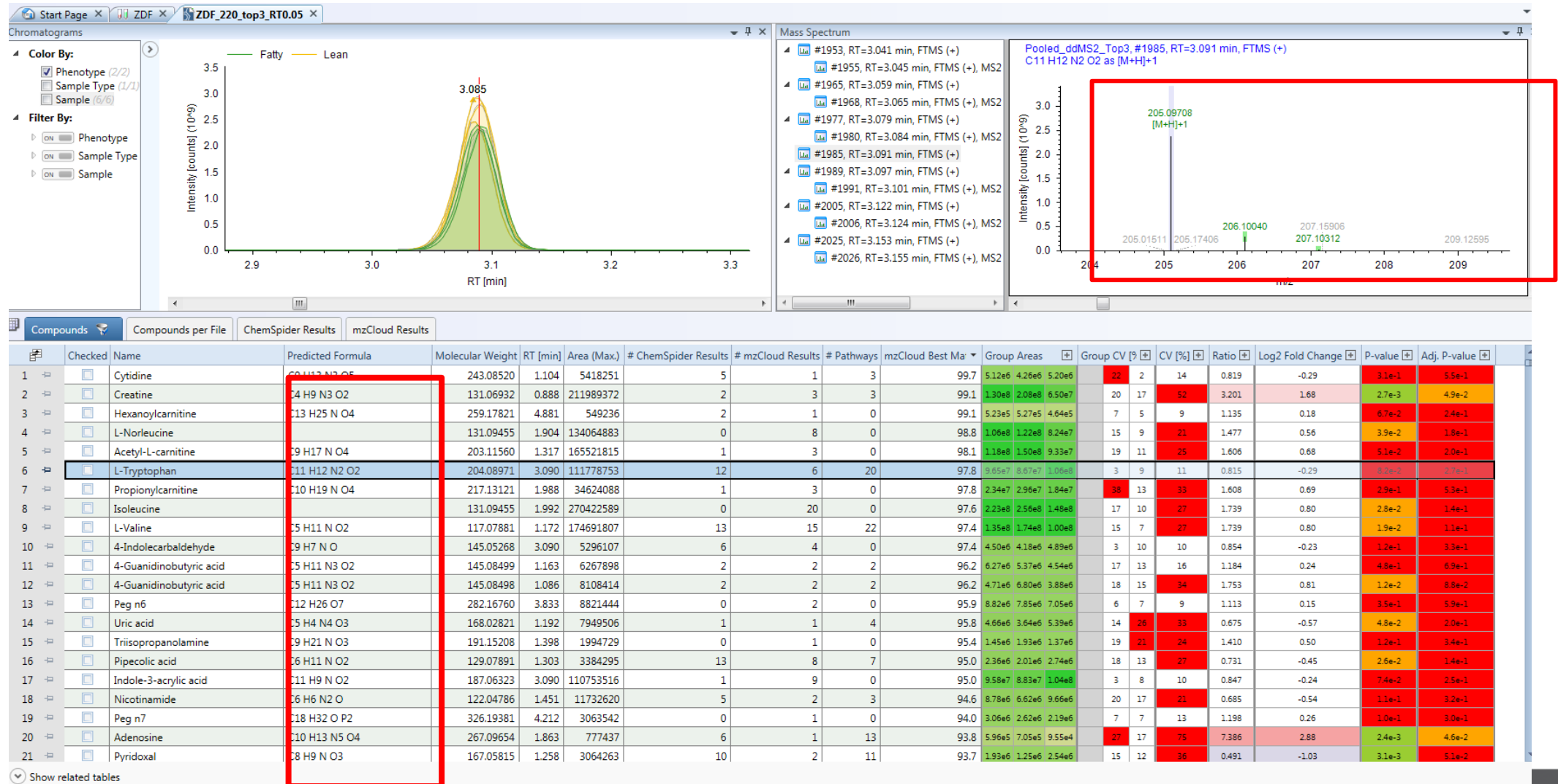

Food Bioactive Compounds



Review

Food Bioactive Compounds and Emerging Techniques for Their Extraction: Polyphenols as a Case Study

José S. Câmara ^{1,2,*}, Bianca R. Albuquerque ^{3,4}, Joselin Aguiar ¹, Rúbia C. G. Corrêa ^{3,5}, João L. Gonçalves ¹, Daniel Granato ⁶, Jorge A. M. Pereira ¹, Lillian Barros ^{3,6} and Isabel C. F. R. Ferreira ³


- ¹ CQM—Centro de Química da Madeira, Campus Universitário da Penteada, Universidade da Madeira, 9020-105 Funchal, Portugal; joselin.aguiar@staff.uma.pt (J.A.); jluis@staff.uma.pt (J.L.G.); jorge.pereira@staff.uma.pt (J.A.M.P.)
- ² Departamento de Química, Faculdade de Ciências Exatas e da Engenharia da Universidade da Madeira, Campus Universitário da Penteada, 9020-105 Funchal, Portugal
- ³ Centro de Investigação de Montanha (CIMO), Instituto Politécnico de Bragança, Campus de Santa Apolónia, 5300-253 Bragança, Portugal; bianca.albuquerque@ipb.pt (B.R.A.); rubiacorreia@ipb.pt (R.C.G.C.); lferreira@ipb.pt (I.C.F.R.F.)
- ⁴ REQUIMTE—Science Chemical Department, Faculty of Pharmacy, University of Porto, Rua Jorge Viterbo Ferreira N° 228, 4050-313 Porto, Portugal
- ⁵ Program of Master in Clean Technologies, Cesumar Institute of Science Technology and Innovation (ICETI), Cesumar University—UnicCesumar, Parana 87050-390, Brazil
- ⁶ Food Processing and Quality, Natural Resources Institute Finland (Luke), Tietotie 2, FI-02150 Espoo, Finland; daniel.granato@luke.fi
- * Correspondence: jsc@staff.uma.pt (J.S.C.); lillian@ipb.pt (L.B.); Tel.: +351-29170-5112 (J.S.C.); +351-2-7333-0901 (L.B.)

Abstract: Experimental studies have provided convincing evidence that food bioactive compounds (FBCs) have a positive biological impact on human health, exerting protective effects against non-communicable diseases (NCD) including cancer and cardiovascular (CVDs), metabolic, and neurodegenerative disorders (NDDs). These benefits have been associated with the presence of secondary metabolites, namely polyphenols, glucosinolates, carotenoids, terpenoids, alkaloids, saponins, vitamins, and fibres, among others, derived from their antioxidant, antiatherogenic, anti-inflammatory, antimicrobial, antithrombotic, cardioprotective, and vasodilator properties. Polyphenols as one of the most abundant classes of bioactive compounds present in plant-based foods emerge as a promising approach for the development of efficacious preventive agents against NCDs with reduced side effects. The aim of this review is to present comprehensive and deep insights into the potential of polyphenols, from their chemical structure classification and biosynthesis to preventive effects on NCDs, namely cancer, CVDs, and NDDs. The challenge of polyphenols bioavailability and bioaccessibility will be explored in addition to useful industrial and environmental applications. Advanced and emerging extraction techniques will be highlighted and the high-resolution analytical techniques used for FBCs characterization, identification, and quantification will be considered.

Keywords: food bioactive compounds (FBCs); polyphenols; disease protection; cancer; cardiovascular diseases (CVDs); neurodegenerative diseases (NDs); bioavailability and bioaccessibility

1. Introduction

Non-communicable diseases (NCDs), mainly cardiovascular diseases (CVDs), cancer, chronic respiratory diseases, neurodegenerative diseases (NDs), and diabetes, represent a severe burden worldwide. According to the most recent data, NCDs caused over 70% of the 57 million deaths worldwide in 2016 [1]. The major risk factors contributing to this scenario have been identified as the combination of an unhealthy diet, poor physical activity, and alcohol and smoking abuse. Diet, in particular, is considered a leading risk factor for illness, death, and disability and it is estimated that one in five deaths are associated with a

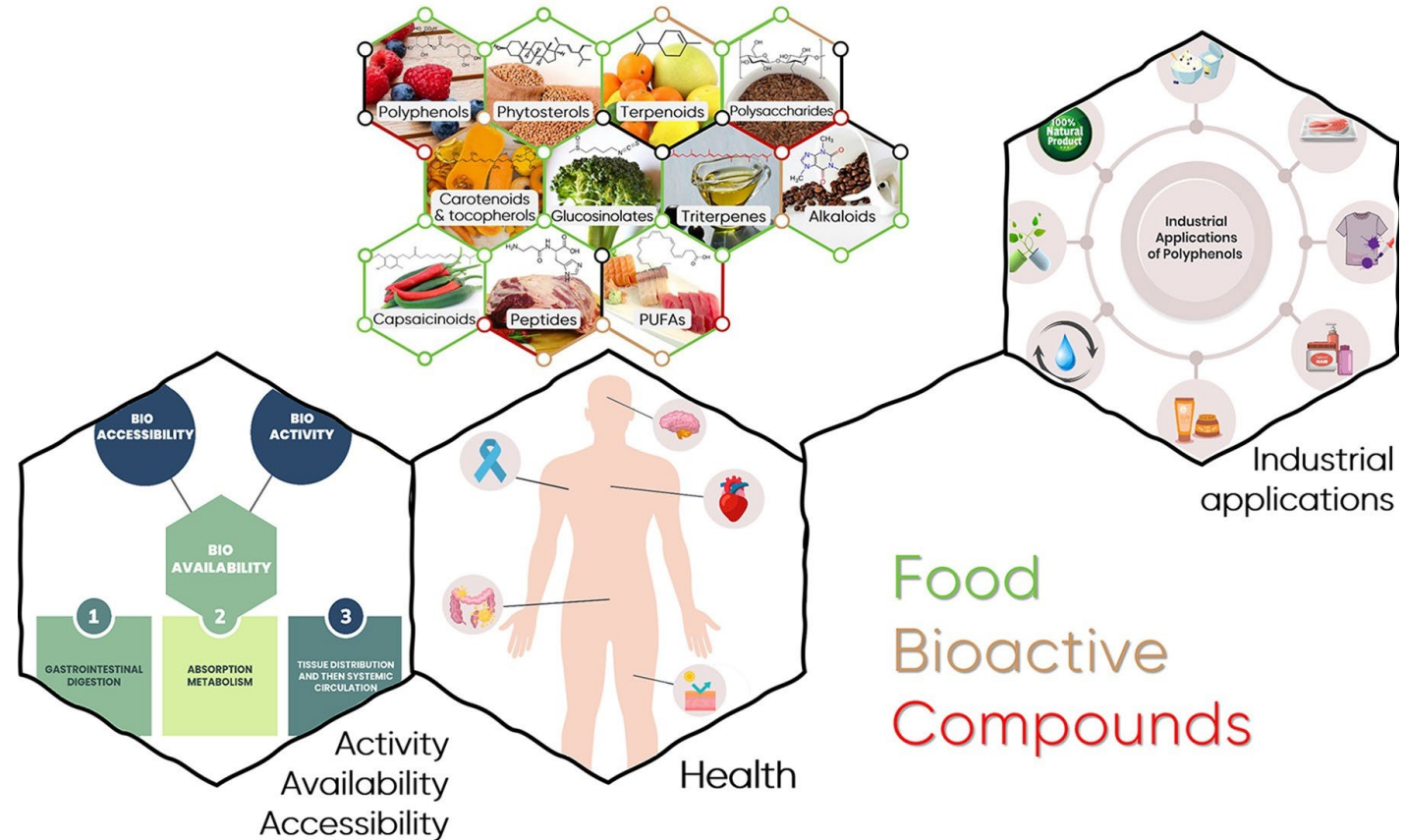

Citation: Câmara, J.S.; Albuquerque, B.R.; Aguiar, J.; Corrêa, R.C.G.; Gonçalves, J.L.; Granato, D.; Pereira, J.A.M.; Barros, L.; Ferreira, I.C.F.R. Food Bioactive Compounds and Emerging Techniques for Their Extraction: Polyphenols as a Case Study. *Foods* **2021**, *10*, 37. <https://doi.org/10.3390/foods1001037>

Received: 27 November 2020
Accepted: 21 December 2020
Published: 24 December 2020

Publisher's Note: MDPI stays neutral with regard to jurisdictional claims in published maps and institutional affiliations.



Copyright: © 2020 by the authors. Licensee MDPI, Basel, Switzerland. This article is an open access article distributed under the terms and conditions of the Creative Commons Attribution (CC BY) license (<https://creativecommons.org/licenses/by/4.0/>).



Orbitrap HRMS in Food Analysis

Food Additives & Contaminants: Part A, 2015
Vol. 32, No. 10, 1568–1606, <http://dx.doi.org/10.1080/19440049.2015.1057240>



Future perspectives in Orbitrap™-high-resolution mass spectrometry in food analysis: a review

Hamide Z. Senyuva^{a*}, Vural Gökmen^b and Ebru Ates Sarikaya^c

^aFoodLife International Ltd, ODTU Teknokent, 06800 Ankara, Turkey; ^bDepartment of Food Engineering, Hacettepe University, 06800 Ankara, Turkey; ^cThermo Fisher Scientific, Dreieich, Germany

(Received 11 March 2015; accepted 28 May 2015)

A literature search from 2007 to 2014 was conducted to identify publications where principally LC-Orbitrap™-high-resolution mass spectrometry (HRMS) has been employed in food analysis. Of a total of 212 relevant references, only 22 papers were from 2007–10, but in subsequent years there has been a steady growth in publications with 38–55 relevant papers being published each year from 2011 to 2014. In the food safety area, over 50% of the published papers were equally divided between pesticides, veterinary drug residues and natural toxins (including mycotoxins) focused primarily on multi-analyte target analysis. LC-Orbitrap-HRMS was also found to be increasingly important for the analysis of bioactive substances, principally phenolic compounds in foods. A number of studies reported for the first time the identification of new fungal metabolites, predominantly various conjugated forms of known mycotoxins. Novel process contaminants were also identified by LC-Orbitrap-HRMS, as were various substances used for food adulteration and bioactive substances in herbal products and dietary supplements. Untargeted analysis is seen as a major future trend where HRMS plays a significant role. Retrospective analysis of scanned high-resolution mass spectra in conjunction with relevant databases can provide new insights. Metabolomics is also being increasingly used where foods are being profiled through fingerprinting using HRMS. All evidence points towards future growth in the number of applications of HRMS in food safety and quality, as the power of this technique gains wider recognition.


Keywords: review; Orbitrap; high resolution mass spectrometry; pesticides; veterinary drugs; natural toxins; process contaminants; adulteration; bioactive substances; non-target analysis; metabolomics

Introduction

Over the past 20 years there have been a number of key developments in mass spectrometry (MS) which have led to step-changes in the way food analysis has been conducted. In LC-MS the development of electrospray (ESI) and atmospheric pressure chemical ionisation (APCI) interfaces together with progressive improvements in MS designs opened up LC-MS to the wider analytical community. Improvements in the ease of use of LC-MS through better software, improved reliability and lower purchase costs have led to significantly wider accessibility, which had been initially restricted to specialists. Lately, further improvements have seen movement from LC-MS to LC-MS/MS, which has now become the standard approach, whether it be with bench-top instruments or more sophisticated MS configurations. Another step change has been the development of high-resolution LC-MS through Orbitrap technology enabling resolutions of 10 000–200 000 to be achieved (Zubarev & Makarov 2013) together with development of time-of-flight (TOF) instruments in the form of LC-TOF-MS in various configurations (Holčápek et al. 2012). The same trend with HRMS, as was the case with LC-MS, has been apparent in the last few years in terms of software being progressively improved and lower capital investment required

with the introduction of LC-HRMS bench-top instruments such as the Exactive™ (a single-stage Orbitrap-MS). There are, what can appear to be, a confusing array of different LC-MS configurations, and a useful overview of commercial systems has been provided comparing the technical specifications of individual manufacturers instruments (Holčápek et al. 2012). Applications of different types of mass detector including Orbitrap™ configurations have been reviewed (Wang, Wang, et al. 2013) highlighting newly developed MS methods, including MALDI TOF-MS (MALDI = matrix-assisted laser desorption/ionisation) imaging and ambient ionisation MS for direct analysis in real time (DART). The advantages and the limitations of different MS techniques in their application to food safety and quality have been assessed (Wang, Wang, et al. 2013).


In the food safety area, the use of HPLC was limited to fluorescence or diode array detection which provided the only viable options offering the sensitivity needed in residue and contaminant analysis. However, with MS available as a detection option, methods quickly migrated onto LC-MS platforms and this transformation was particularly apparent in the pesticide and veterinary drug residue areas, where multi-residue and multi-class methods were developed (Kaufmann et al. 2011a; Mol et al. 2012). A review of LC-MS in food safety

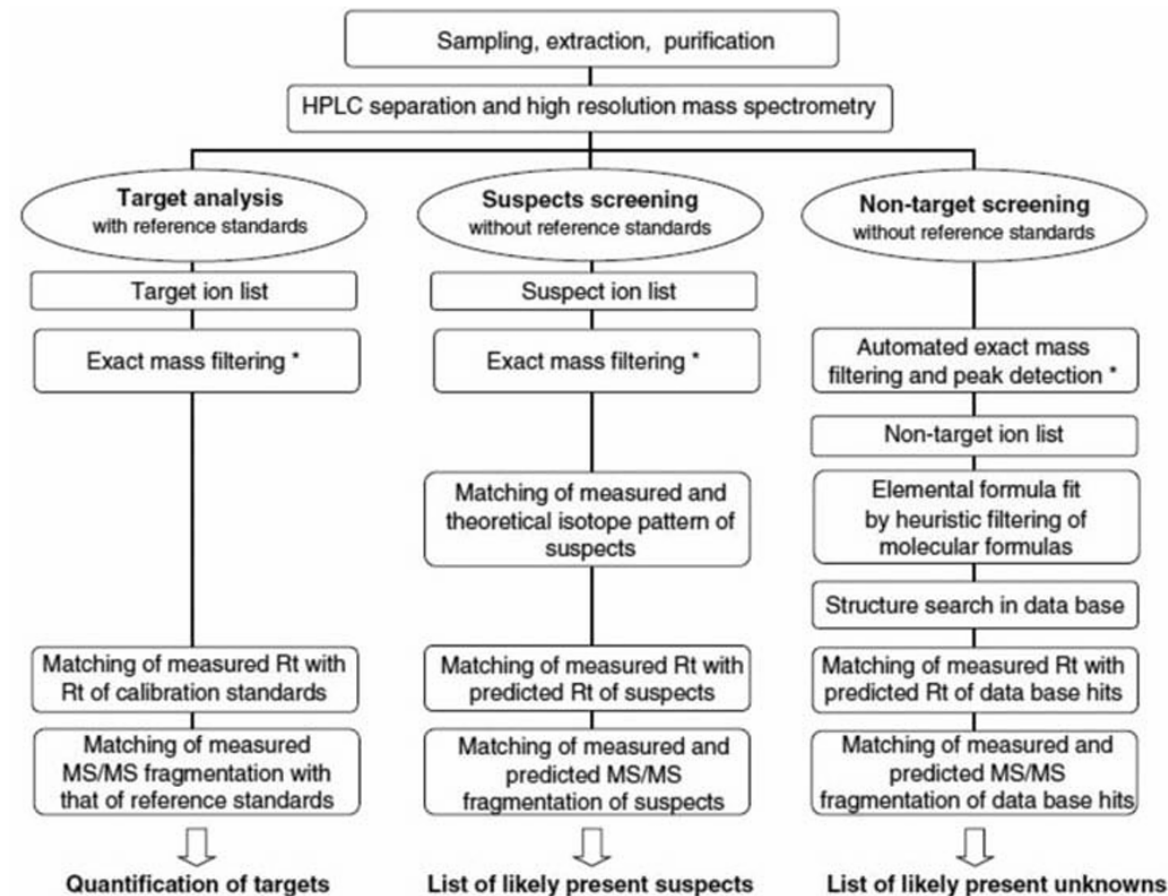

*Corresponding author. Email: hamide.senyuva@foodlifeint.com

Compound Identification

Know more unknowns

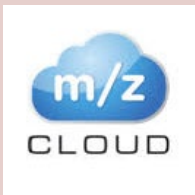Database	Molecular Ion (Precursor Ion)	MS ² Spectrum	Metabolic Pathway
Online	ChemSpider 	mzCloud 	- Metabolika Pathways - KEGG - BioCyc
Offline In-house Database	Mass Lists Optional Import from ChemSpider	mzVault 	Metabolika Pathways

Automatically identify components

- **Spectral library searching**
 - Online mzCloudTM spectral library
 - In-house Thermo ScientificTM mzVaultTM spectral libraries
- **mzCloud can provide similarity search results**, giving you even more information to use for unknown identification
- **Determine elemental composition**, including fine isotopes, using HRAM MS and MS/MS data. Then, use this information to search multiple chemical databases for further annotation information
- **Structure proposals automatically annotated on an experimental spectrum**

Orbitrap HRMS in Food Analysis



Group	Specific compounds (number identified)	Matrix	Basis for identification	Reference
Mycotoxins	Fusarins (10)	Culture medium	LC-HRMS and ¹³ C-NMR	Kleigrew et al. (2012)
	DON conjugates (8)	Wheat	LC-HRMS, SIL	Kluger et al. (2013)
	DON oligoglycosides (3)	Malt, beer, bread	IAC, LC-HRMS	Zachariasova et al. (2012)
	Fusarenon X-glucoside	Wheat	LC-HRMS	Nakagawa et al. (2011)
Marine biotoxins	<i>A. flavus</i> anthraquinone metabolites	Culture medium	LC-HRMS, LC-ITMS	Malysheva et al. (2014)
Plant toxins	Pinnatoxins (6)	Mussels	LC-HRMS and synthesis	McCarron et al. (2012)
	Acetogenins (80)	Tropical fruit	HPLC-UV-LTQ-Orbitrap	Le Ven et al. (2014)
	Quinoline, carboline and glycinamide compounds	Cow's milk	HPLC/DAD/HRMS and 1D and 2D NMR	Rouge et al. (2013)
Veterinary drugs	Enrofloxacin metabolites	Chicken	LC-LTQ-Orbitrap and LC-QqToF	Morales-Gutiérrez et al. (2014)
	Mequindox metabolites (24)	Chicken bile, plasma, faeces, tissues	LC-HRMS	Shan et al. (2012)
Process contaminants	Estradiol-dipeptide biomarker	Urine	LC-LTQ-Orbitrap	Regal et al. (2013)
	MCPD derivatives	Palm oil	LC-LTQ-Orbitrap – mass defect filtering	Nagy et al. (2011)
Adulteration	Mepiquat	Coffee	LC-Q-Exactive MS	Hammel et al. (2014)
	Sildenafil analogue	Dietary supplement	LC-LTQ-Orbitrap-MS	Kee et al. (2013)
Bioactive substances	Basic red 46 dye	Spices	LC-Exactive	Ruf et al. (2012)
	Flavonoids, phenolic acids, phenylpropanoid glycosides (56)	TCM	UHPLC-ESI/LTQ-Orbitrap-MS	Chen et al. (2014)
	Caffeoyl derivatives, flavonoids, coumarins and sesquiterpenoids (50)	TCM	LC-HRMS and ¹ H-NMR, ¹³ C-NMR	Li et al. (2011)
	Flavanol, flavanone, dihydrochalcone derivatives (38 compounds)	Tomatoes	LC-LTQ-Orbitrap-MS and LC-MS/MS	Vallverdu-Queralt et al. (2010)
	Quercotriterpenoside (2 compounds)	Wine	LC-HRMS, LC-MS/MS 2D ¹ H-NMR, ¹³ C-NMR	Marchal et al. (2011)

Food Allergens

Analytical and Bioanalytical Chemistry
<https://doi.org/10.1007/s00216-023-04894-2>

RESEARCH PAPER



In-house validation of an LC–MS method for the multiplexed quantitative determination of total allergenic food in chocolate

Rosa Pilolli¹ · Antonella Lamonaca^{1,2} · Chiara Nitride^{3,4} · Elisabetta De Angelis¹ · Christof van Poucke⁵ · Nathalie Gillard⁶ · Anne-Catherine Huet⁶ · Marc De Loose⁵ · Jean Henrotin⁶ · E. C. N. Mills⁴ · Linda Monaci¹

Received: 18 April 2023 / Revised: 28 July 2023 / Accepted: 31 July 2023
 © The Author(s) 2023

Abstract

Mass spectrometry has been widely accepted as a confirmatory tool for the sensitive detection of undeclared presence of allergenic ingredients. Multiple methods have been developed so far, achieving different levels of sensitivity and robustness, still lacking harmonization of the analytical validation and impairing comparability of results. In this investigation, a quantitative method has been validated in-house for the determination of six allergenic ingredients (cow's milk, hen's egg, peanut, soybean, hazelnut, and almond) in a chocolate-based matrix. The latter has been produced in a food pilot plant to provide a real and well-characterized matrix for proper assessment of method performance characteristics according to official guidelines. In particular, recent considerations issued by the European Committee for Standardization have been followed to guide a rigorous single-laboratory validation and to feature the main method performance, such as selectivity, linearity, and sensitivity. Synthetic surrogates of the peptide markers have been used both in native and labelled forms in matrix-matched calibration curves as external calibrants and internal standards, respectively. A two-order of magnitude range was investigated, focusing on the low concentration range for proper assessment of the detection and quantification limits (LOD and LOQ) by rigorous calibration approach. Conversion factors for all six allergenic ingredients have been determined for the first time to report the final quantitative information as fraction of total allergenic food protein (TAFP) per mass of food ($\mu\text{g}_{\text{TAFP}}/\text{g}_{\text{food}}$), since such a reporting unit is exploitable in allergenic risk assessment plans. The method achieved good sensitivity with LOD values ranging between 0.08 and 0.2 $\mu\text{g}_{\text{TAFP}}/\text{g}_{\text{food}}$ for all ingredients besides egg and soybean, whose quantitative markers reported a slightly higher limit (1.1 and 1.2 $\mu\text{g}_{\text{TAFP}}/\text{g}_{\text{food}}$, respectively). Different samples of chocolate bar incurred at four defined concentration levels close to the currently available threshold doses have been analyzed to test the quantitative performance of the analytical method, with a proper estimate of the measurement uncertainty from different sources of variability. The sensitivity achieved resulted in compliance with the various threshold doses issued or recommended worldwide.

Keywords Food allergen · Mass spectrometry · Quantitative method · In-house validation · Uncertainty · Conversion factors

Highlights of Orbitrap Mass Spectrometry in Allergen Detection

-  Simultaneous detection of multiple allergens
 → Covers 5 types: low-fat milk, whole egg, soy flour, hazelnut, and peanut.
-  Protein extraction via ultrasonic treatment
 → Simple 30-minute protocol + cleanup using exclusion cartridge.
-  Full MS/DIA acquisition mode
 → Ensures high sensitivity and precision in measurement.
-  17 allergen-specific peptides detected
 → All identified in a single run from processed cookie samples.
-  Detection limits: 60–100 $\mu\text{g}/\text{g}$
 → Capable of detecting trace levels in complex food matrices.
-   Validated with in-lab precision testing
 → Demonstrates reliability and consistency for real-world use.
-  Superior to traditional methods (ELISA / PCR)
 → No need for allergen-specific kits
 → Saves time, expands coverage
 → Detects processed allergens with altered protein structures

Protein & Peptide Biomarkers



Available online at www.sciencedirect.com

ScienceDirect

Current Opinion in
Food
Science

Proteins and peptides: proteomics approaches for food authentication and allergen profiling

Mónica Carrera¹, Ana G Abril¹, Manuel Pazos¹, Pilar Calo-Mata², Tomás G Villa³ and Jorge Barros-Velázquez²

Food authentication and food allergy profiling are relevant topics that must be carefully considered by authorities, food production and manufacturing industries, and customers. Precise, sensitive, and efficient detection procedures are crucial for authenticating food and identifying allergens. In this paper, we present a current review of advanced proteomic approaches for food authentication and allergen profiling. We discuss three proteomics strategies, namely, discovery proteomics, targeted proteomics, and proteomics-based systems biology, used to authenticate food and profile allergens. This review includes a selection of the most representative papers published in this field in the last few years. Furthermore, we discuss future directions and potential opportunities in this field.

Addresses

¹Spanish National Research Council (CSIC), Institute of Marine Research (IM-CSIC), Department of Food Technology, 36208 Vigo, Spain

²School of Veterinary Sciences, University of Santiago de Compostela, Campus Lugo, Department of Analytical Chemistry, Nutrition and Food Science, Food Technology Division, 27002 Lugo, Spain

³Faculty of Pharmacy, University of Santiago de Compostela, Department of Microbiology and Parasitology, 15898 Santiago de Compostela, Spain

Corresponding author: Carrera, Mónica (mcarrera@im.csic.es)

Current Opinion in Food Science 2024, 57:101172

This review comes from a themed issue on **Foodomics Technologies**

Edited by Markus Fischer

For complete overview of the section, please refer to the article collection, "Foodomics Technologies 2024"
Available online 6 May 2024

<https://doi.org/10.1016/j.cofs.2024.101172>

2214-7993/© 2024 The Author(s). Published by Elsevier Ltd. This is an open access article under the CC BY-NC-ND license (<http://creativecommons.org/licenses/by-nc-nd/4.0/>).

Introduction


Food authentication and allergy profiling are important measures to ensure food quality and safety. Confirming the authenticity of food products is essential for preventing the consumption of harmful substances and

protecting consumer health. Regulatory bodies are increasingly focused on accurately verifying products and preventing the substitution of species, addressing the concerns of customers and the food industry. These regulations comprise the U.S. Food and Drug Administration's Food Safety Modernization Act 2011, the European Regulation (EC) No 178/2002 identified as the General Food Law Regulation, and the Food Information to Consumers Regulation (Regulation [EU] No 1169/2011). The guidelines require food labels to clearly state the species or commercial name of the product.

Accurate labels with information about potential allergens also help people with food allergies obtain safe food. Approximately 6–8% of kids and 2–4% of adults are diagnosed with food allergies [1]. The European Food Safety Authority has recognized 14 food items as allergens, including fish, mollusks, shellfish, milk, eggs, soybeans, nuts, celery, peanuts, mustard, wheat, lupin, sesame, and sulfur dioxide/sulfites. Numerous countries have implemented stringent regulations for the labeling of allergens to safeguard consumers, as outlined in Directive 2007/68/EC.

Authorities, the food industry, and customers play a role in guaranteeing food quality/safety, as food authentication and allergy profiling have become topics of interest. To meet these guidelines, accurate and efficient detection methods are needed for identifying food species and controlling allergies. Various analytical methodologies have been applied to authenticate food and profile allergens. There are several commonly utilized methods, including high-performance liquid chromatography, near-infrared spectroscopy, nuclear magnetic resonance spectroscopy, immunoassays, and stable isotope ratio; in addition, genomics, metabolomics, and proteomics approaches have been recently applied [2–4]. In this sense, omics methodologies have profoundly impacted the research community and are increasingly being utilized in modern food science.

Research in the proteomics field examines proteins within a biological system on a large scale [5]. The objective of this field is to clarify the assembly, function, modifications, interactions, and regulation of proteins, thereby offering valuable knowledge on the dynamic

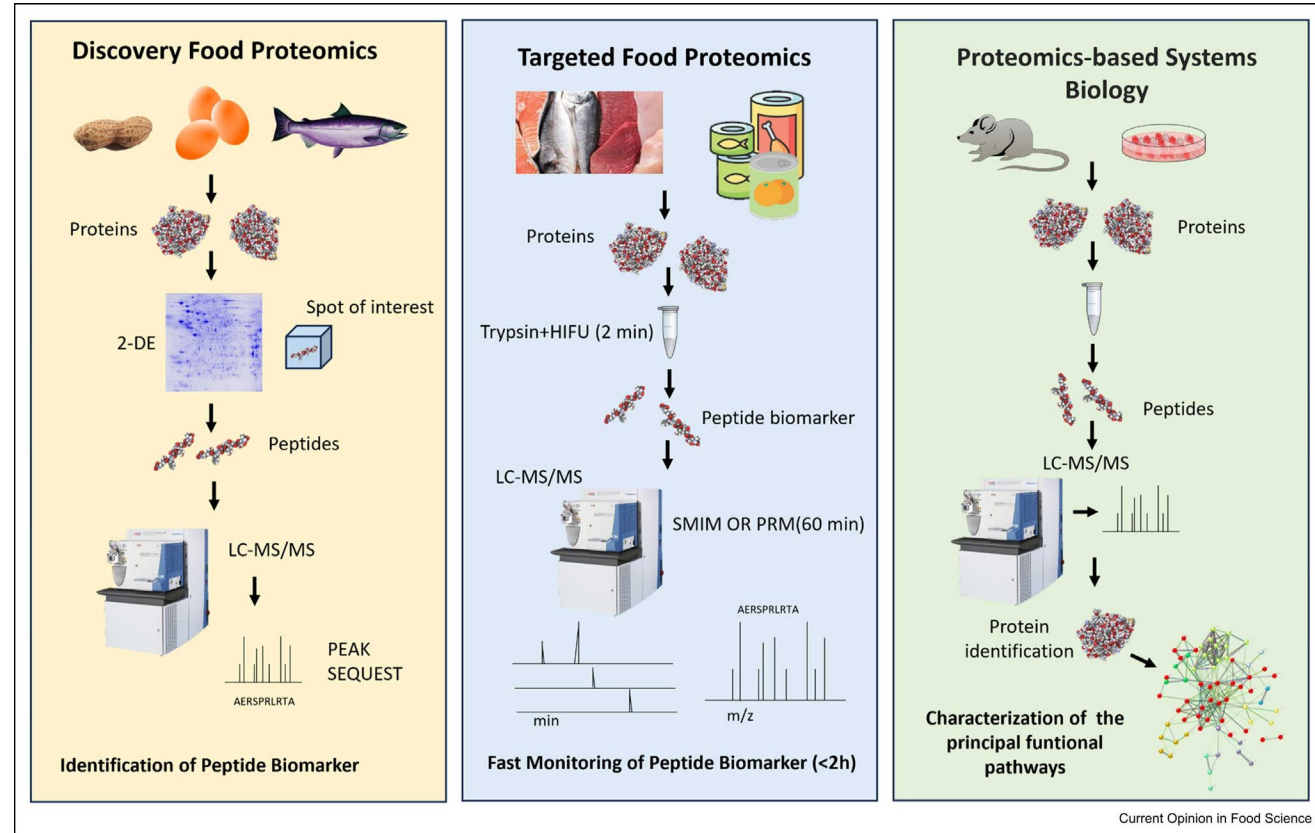

Current Opinion in Food Science